


Quad Tc/mV Amp

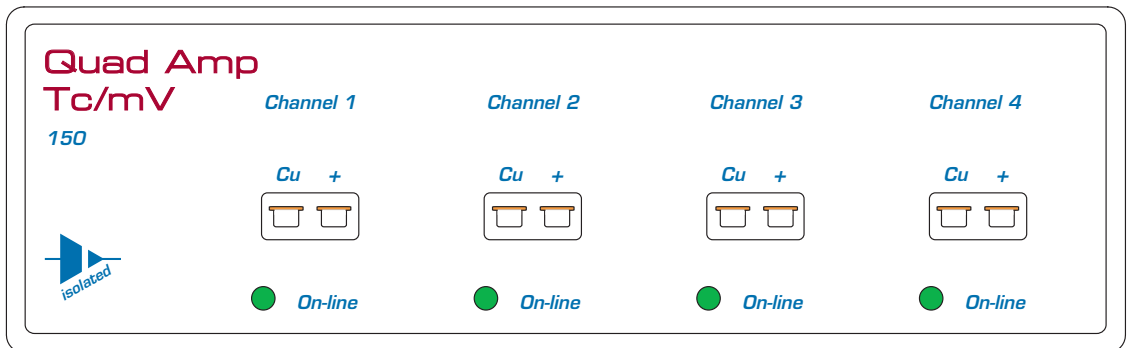

This document was, as far as possible, accurate at the time of printing. Changes may have been made to the software and hardware it describes since then: eDAQ Pty Ltd reserves the right to alter specifications as required. Late-breaking information may be supplied separately. Latest information and software updates can be obtained from our web site.

Trademarks of eDAQ

e-corder and PowerChrom are registered trademarks of eDAQ Pty Ltd. Specific model names of data recording units, such as e-corder 201, and PowerChrom 280, are trademarks of eDAQ Pty Ltd. Chart and Scope are trademarks of ADInstruments Pty Ltd and are used under license by eDAQ. EChem is a trademark of eDAQ Pty Ltd.

Other Trademarks

Mac OS, and Macintosh, are registered trademarks of Apple Computer, Inc. Windows 98, Windows Me, Windows 2000, and Windows XP are trademarks of Microsoft Corporation.

PostScript, and Acrobat are registered trademarks of Adobe Systems, Incorporated.

All other trademarks are the properties of their respective owners.

Product: Quad Tc/mV Amp (EA150)

Document Number: U-EA150-0806


Copyright © August 2006

eDAQ Pty Ltd
6 Doig Avenue
Denistone East, NSW 2112
AUSTRALIA

<http://www.eDAQ.com>
email: info@eDAQ.com

All rights reserved. No part of this document may be reproduced by any means without the prior written permission of eDAQ Pty Ltd.



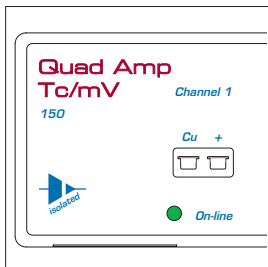

Contents

1 Overview 1

How to Use this Manual 2
Checking the Quad Tc/mV Amp 2
eDAQ Amps 2

2 The Quad Tc/mV Amp 5

The Front Panel 6
 The Thermocouple Inputs 6
 The On-line Indicators 6
The Back Panel 7
 The Signal Outputs 7
 I²C Connector 7
Connecting to the e-corder 7
 The Quad Tc/mV Amp Self Test 9
Using Chart software 9
 Setting up the Quad Tc/mV Amp 9
 Previewing the Signal 10
Calibration 12
 Automatic Calibration 12
 13
 Manual Calibration 13

A Technical Aspects 15

Quad Tc/mV Amp Operation 15
Technical Description 15

B Troubleshooting 17

C Specifications 19

Index 21

License & Warranty 23

1

CHAPTER ONE

Overview

The eDAQ Quad Tc/mV Amp is designed to work with an e-corder system, recording signals from up to four thermocouples (or other low voltage sources) simultaneously.

How to Use this Manual

This manual describes how to set up and begin using your Quad Tc/mV Amp. The use of the pH/mV Amp with the Chart software Thermocouple Calibration extension is discussed. The appendices provide technical and troubleshooting information.

Checking the Quad Tc/mV Amp

Before you begin working with the Quad Tc/mV Amp, please check that:

- all items described in the packing list are included; and that
- there are no signs of damage that may have occurred during transit.

Contact your eDAQ distributor if you encounter a problem.

You should also become familiar with the basic features of your e-corder system, which are discussed in the *e-corder Manual* which will be installed on your hard disk.

eDAQ Amps

The Quad Tc/mV Amp is part of a family of preamplifiers known as eDAQ Amps.

The Quad Tc/mV Amp is designed for recording signals from up to four pH and other potentiometric electrodes. As with other eDAQ Amps, it is designed to be operated under full software control and is automatically recognized by Chart or Scope software which controls its gain range, signal filtering, and other settings.

The range of eDAQ Amps includes the:

- **Quad pH/mV Amp**, suitable for connection of pH, ion selective, and potentiometric (ORP) electrodes
- **Potentiostat**, a three-electrode potentiostat that can be used for voltammetric and amperometric experiments. Gain ranges of 20 nA to 100 mA in 1:2:5 steps.
- **Picostat**, a high sensitivity three-electrode potentiostat suitable for use with carbon fibre and other microelectrodes. Current gain ranges of 10 pA to 100 nA in 1:2:5 steps.
- **QuadStat**, a four channel, three electrode potentiostat suitable for use with amperometric biosensors. Current gain ranges of 2 nA to 1 mA/channel in 1:2:5 steps.
- **Bridge Amp**, suitable for sensors requiring a low drift, high gain differential amplifier. Also provides DC excitation
- **GP Amp**, suitable for high output sensors requiring a high impedance single ended or differential amplifier. Also provides DC excitation.

See our web site at www.eDAQ.com for more information.



2

C H A P T E R T W O

The Quad Tc/mV Amp

This chapter describes how to connect and use your eDAQ Quad Tc/mV Amp.

IMPORTANT: Always make sure that the e-corder is turned off before you connect or disconnect the Quad Tc/mV Amp. Failure to do this may result in damage to the e-corder and/or the Quad Tc/mV Amp.

The Front Panel

The front panel of the Quad Tc/mV Amp is shown in Figure 2-1.

The Thermocouple Inputs

The miniature thermocouple input connectors marked 'Cu' and '+' of the Quad Tc/mV Amp provide connection for most types of thermocouple.

The On-line Indicators

When lit, these indicate that the Chart software has located and initialized the Quad Tc/mV Amp. If the light does not go on when the software is run, check that the Quad Tc/mV Amp is properly connected.

Figure 2-1
The Quad Tc/mV Amp front panel


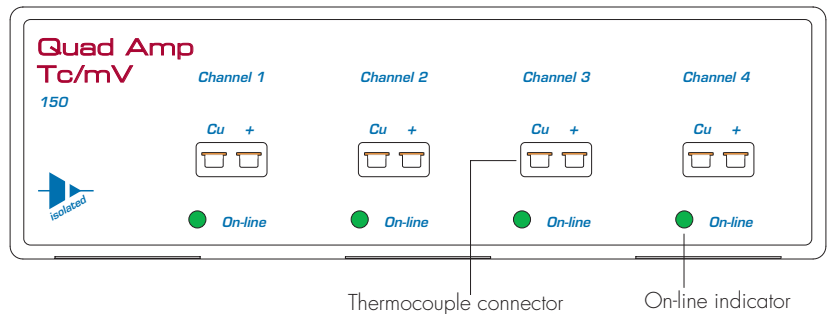


Figure 2-2
The Quad Tc/mV Amp back panel

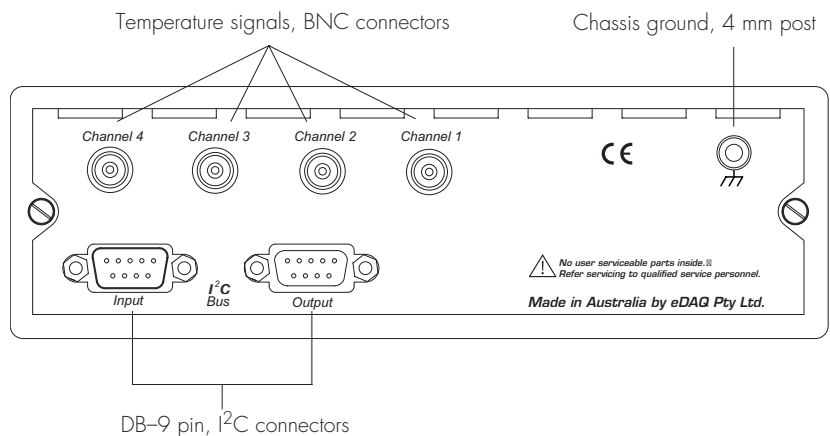

The Back Panel

The back panel of the Quad Tc/mV Amp is shown in [Figure 2-2](#).

The Signal Outputs

The Quad Tc/mV Amp provides four signals, one for each sensing thermocouple. These signals are accessed on the Quad Tc/mV Amp back panel via four BNC connectors labelled Channel 1 — Channel 4.


I²C Connector

The back panel, [Figure 2-2](#), has two DB-9 pin 'I²C bus' connectors labelled Input and Output. The Input connector enables connection to the e-corder (or to the output of other eDAQ Amps). This connection provides power to the Quad Tc/mV Amp and carries the various control signals (for gain range and filter selection) to and from the e-corder. The pin assignments are shown in [Figure 2-3](#).

Connecting to the e-corder

Your Quad Tc/mV Amp will have been supplied with an I²C cable (DB-9 pin connectors at either end), and four cables with BNC connectors at either end.

Figure 2-3
I²C connector pin assignments

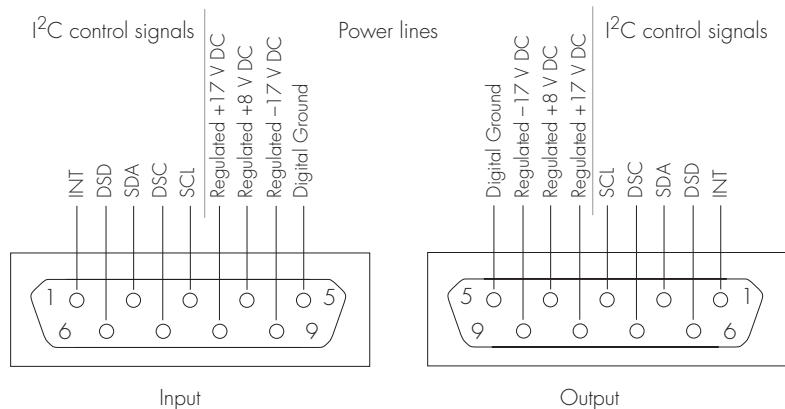

First make sure that the e-corder is turned off. Then connect the I²C cable to the I²C connector on the back panel of the e-corder, and the other end to the I²C Input connector on the back panel of the pH/mV Amp. Use BNC cables to connect the signal outputs on the back panel of the Quad Tc/mV Amp to the front panel of the e-corder as shown in Figure 2–4 and Figure 2–5.

Check that all connections are firm. Loose connections may cause failure to recognize the eDAQ Amp by the software, erratic behavior, or loss of signal.

Multiple eDAQ Amps can be connected to an e-corder. The number that can be connected depends on the number of input channels on the e-corder. The initial eDAQ Amp should be connected as shown in Figure 2–5. The remainder are linked via I²C cables, connecting the I²C output of one eDAQ Amp to the I²C input of the next. The signal outputs of each eDAQ Amp are connected to an input channel of the e-corder.

Figure 2–4

Four thermocouples shown connected to a Quad Tc/mV Amp and e-corder 401


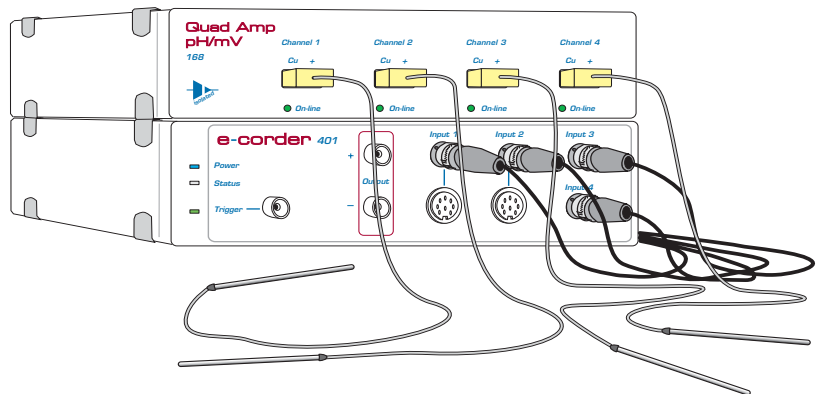


Figure 2–5

Back view of Quad Tc/mV Amp connected to e-corder

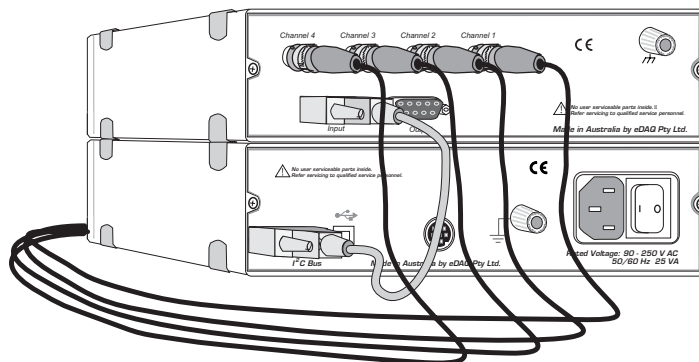

The Quad Tc/mV Amp Self Test

After you have installed the software, connected the e-corder to a computer as described in your *e-corder Manual*, and connected the Quad Tc/mV Amp as described above, a quick check can be performed:

- Turn on the e-corder and check that it performs its normal diagnostic tests, as described in the *e-corder Manual*
- Once the e-corder is ready, open Chart software.
- During initialization, you should see these indicators flash briefly and then remain lit — the Quad Tc/mV Amp is ready for use.


If the indicators do not light, check your cable connections and repeat the procedure. If this does not solve the problem, contact your eDAQ distributor.

Using Chart software

Setting up the Quad Tc/mV Amp

When the Quad pH/mV Amp is connected to an e-corder input and successfully installed, the Quad Thermocouple... menu command appears in the Chart software Channel Function pop-up menu, [Figure 2-6](#).

Figure 2-6
Accessing the Quad Tc/mV Amp controls

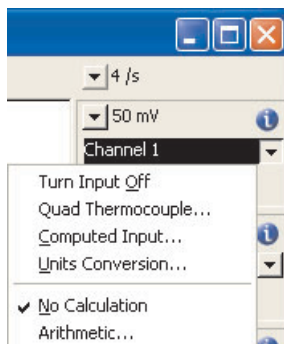


Choosing the Quad Thermocouple... menu command will access the Quad Thermocouple dialog box (which replaces the normal e-corder

Input Amplifier dialog box for that e-corder channel — your *Chart Software Manual* have details on the Input Amplifier dialog box).

Previewing the Signal

The Quad Thermocouple dialog box, [Figure 2–6](#), allows you to preview a signal so that you can select the amplification and filter settings.

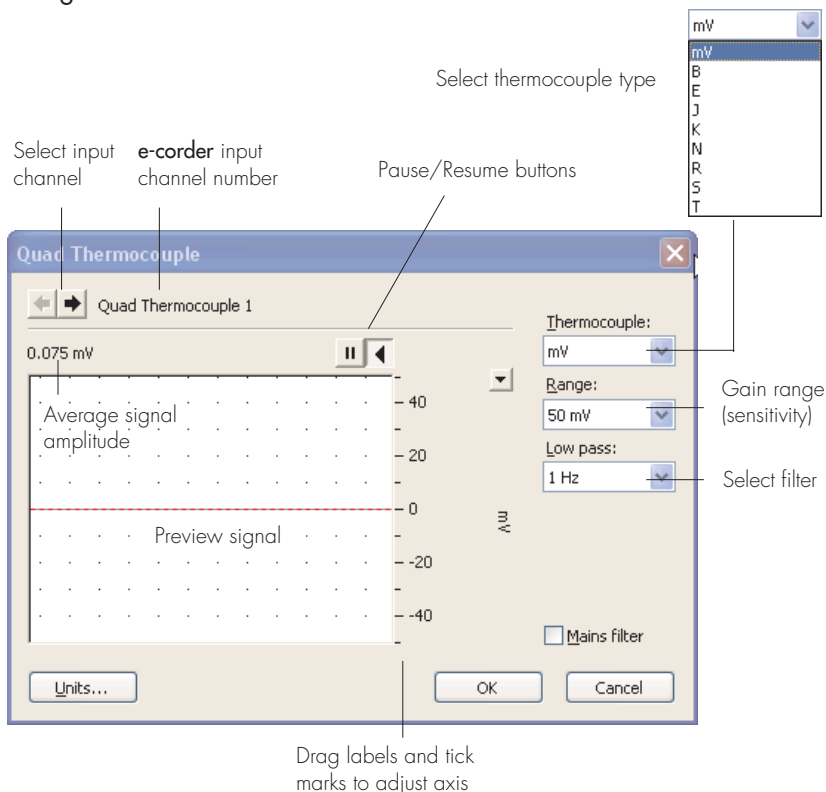

Figure 2–7
The Quad Tc/mV Amp dialog box, for previewing a signal (Chart software on a Windows Computer)



The incoming signal is displayed in real time, but is not recorded to hard disk (once the signal moves across the display area it is lost). Click the OK button to apply the selected settings.

Signal Display

The input signal is displayed so that you can see the effect of changing the settings. The average signal value is displayed at the top left of the display area.

You can stop the signal scrolling by clicking the Pause button at the top right of the display area . Click the Scroll button  to resume scrolling.

You can shift and stretch the vertical Amplitude axis to make the best use of the available display area (see the Input Amplifier dialog in the *Chart Software Manual* for further details). Changes made here update settings in the main window of the program.

Thermocouple Type

Select the appropriate thermocouple type from this menu. If 'mV' is chosen then the raw millivolt signal will be shown. If any other selection is made then the ice point corrected thermocouple signal will be shown in units of mV corresponding to the emf values listed in the NIST ITS-90 thermocouple tables, see <http://srdata.nist.gov/its90/main/>.

Setting the Range

Use the Range pop-up menu to select the input range (sensitivity) of the channel. Available ranges are shown in [Table 2-1](#). Signals in excess of the range value will be truncated.

Filtering


The low pass filter will remove high frequency noise. For most thermocouple probes filter values of 1 – 10 Hz will be appropriate.

The Mains Filter, Mains filter, removes 50 or 60 Hz interference due to mains hum. See the *Chart Software Manual* for more information. For most applications it is recommended to have the Mains Filter on.

Table 2-1
Ranges available for different input settings.

mV	Available ranges (°C) for various thermocouple types							
	B	E	J	K	N	R	S	T
±20	1820	290	360	485	585	1680	1760	250
±10	1500	150	180	250	320	960	1040	115
±5	1020	80	95	120	170	550	580	60
±2	630	35	40	50	75	260	265	30

Units Calibration

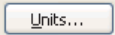Clicking the  button access the Units Conversion dialog box. For more information about Units Conversion, see the *Chart Software Manual*. This type of calibration should only be used for small temperature ranges that can be calibrated with only two accurately known temperature sources.

See the next section for calibration of probes, without the need of accurately known temperature sources.

When calibration is required over a wide range of temperature and several accurately known temperature sources are to be employed then use the Multiple Point Calibration extension.

Calibration

Automatic Calibration

The Thermocouple extension is a Chart software plug-in that adds the 'Thermocouple...' command to the Chart channel pop-up menus. The latest version of the extension is available at http://www.edaq.com/chart_extns_list.html

This extension:

- allows the user to select the type of thermocouple (B, E, J, N, K, R, S or T) being used;
- applies a polynomial correction to convert the voltage signal to degrees (°C, °F, or K) according to the thermocouple ITS-90 tables, see <http://srdata.nist.gov/its90/main/>

The resulting temperature values will usually be accurate within the tolerance of the thermocouple being used — consult your thermocouple probe documentation. In most cases thermocouples conform to either the ANST/ANSI E230 Standard, or Special, Limits of Error, see [Table 2-2](#). In some cases, especially over limited temperature ranges, manufacturers may guarantee better tolerances.

Table 2-2
Limits of Error for
Thermocouples,
Reference Junction
0 °C (32 °F)

Tc Type	Temperature Range		Limits of Error	
	°C	°F	Standard	Special
T	-200 to -66	-328 to -87	±1.5%	—
	-66 to 0	-87 to +32	±1°C, ±2°F	—
	0 – 133	32 – 270	±1°C, ±2°F	±0.5°C, ±1°F
	133 – 350	270 – 662	±0.75%	±0.4%
J	0 – 293	32 – 559	±2.2°C, ±4°F	±1.1°C, ±2°F
	293 – 750	559 – 1382	±0.75	0.4%
E	-200 to -100	-328 to -148	±1%	—
	-100 to 0	-148 to +32	±1.1°C, ±3°C	—
	0 – 340	32 – 644	±1.7°C, ±3°F	±1°C, ±2°F
	340 – 900	644 – 1652	±0.5%	±0.4%
K	-200 to -110	-328 to -166	±2%	—
	-110 to 0	-166 to +32	±2.2°C, ±4°F	—
	0 – 293	32 – 559	±2.2°C, ±4°F	±1.1°C, ±2°F
	293 – 1250	559 – 2282	±0.75%	±0.4%
N	0 – 293	32 – 559	±2.2°C, ±4°F	±1.1°C, ±2°F
	0 – 1250	559 – 2282	±0.75%	±0.4%
R, S	0 – 600	32 – 1112	±1.5°C, ±3°C	±0.6°C, ±1°F
	600 – 1450	1112 – 2642	±0.25%	±0.1%
B	870 – 1700	1598 – 3092	±0.5%	—

Manual Calibration

If known temperature sources are available then very accurate calibration can be achieved using the raw mV signal setting. The signal can be calibrated over small ranges with a linear function using the Chart software Units Conversion feature (see the *Chart Software Manual*). Over wider ranges the Multipoint Calibration extension software plug-in can be used, see http://www.edaq.com/chart_extns_list.html



A

A P P E N D I X A

Technical Aspects

This appendix describes some of the important technical aspects of the Quad Tc/mV Amp. You do not need to know the material here to use the Tc/mV Amp. Please remember that user modification of the equipment voids your rights under warranty.

Quad Tc/mV Amp Operation

The Quad Tc/mV Amp, and other eDAQ Amps, have been designed to be used with an e-corder system. The Tc/mV Amp provides:

- high impedance, electrically isolated inputs suitable for direct connection to most thermocouples;
- cold junction compensation;
- software controlled filtering of the signal;
- software selectable gain ranges.


Technical Description

As with other eDAQ Amps, all internal functions of the Quad Tc/mV Amp are controlled from the e-corder by sending information on a special communications connection called the I²C (eye-squared-sea) bus. This connection supplies power to the eDAQ Amp as well. eDAQ Amps are connected to the analog input channels of the e-corder via BNC-to-BNC cables, through which the amplified and filtered signal is

sent to the e-corder. The overall operation of the pH/mV Amp can be better understood by referring to [Figure A-1](#).

The thermocouple signal is first amplified ($\times 100$) and then passed through an isolation transformer. The signal is then ice point corrected, low pass filtered and then sent to the e-corder for secondary amplification and filtering.

Figure A-1
Block Diagram of the
Quad Tc/mV Amp

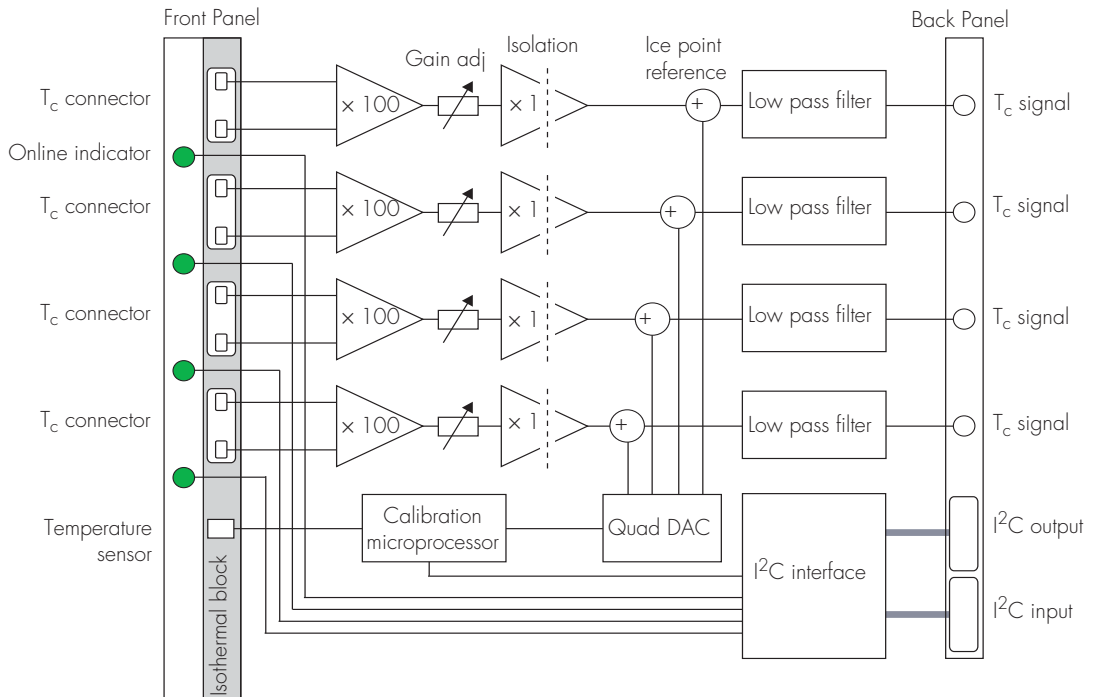

B

A P P E N D I X B

Troubleshooting

If you experience any difficulty with the Quad pH/mV Amp use this section to try to resolve the problem. In the majority of cases, the problem can usually be fixed by ensuring that the connecting cables are correctly connected and firmly attached and the re starting the Chart or Scope software.

The On-Line indicator fails to light when Chart or Scope is opened

- If the e-corder is off check switches and power connections.
- The BNC to BNC cables from the Tc/mV Amp to the analog input channels of the e-corder may not be properly connected. Check that each cable is firmly connected to the BNC connector on the back panel of the pH/mV Amp and to a BNC input on the front panel of the e-corder.

All or some of on-line indicators fail to light when the Chart software is opened, and the pH/mV Amp controls are not accessible

- If all the lights are off, then the I²C cable from the Tc/mV Amp to the e-corder may not be properly connected. Check to see that the I²C cables are firmly seated and screwed in.
- Try a new cable to check that the BNC and/or I²C cables are working correctly.
- If you have tried the above suggestions and still cannot get the Quad Tc/mV Amp to work properly, then try using it on another e-corder if you have access to one. If the on-line indicators still fail to light on the second e-corder, the Quad Tc/mV Amp unit may be faulty and should be returned for repair.



C

A P P E N D I X C

Specifications

Quad Tc/mV Amp

Tc/mV Amplifier:	Four. High impedance, electrometer type. Galvanically isolated.
Input range:	± 20 mV
Secondary ranges:	$\pm 10, 5, 2$ mV, supplied by e-corder
Input resolution:	16 bits, 0.0015% of range
Input impedance:	$10^{13} \Omega$ typical
Amplifier noise:	$< 1 \mu\text{V rms}$ ($4 \mu\text{V p-p}$) @ DC – 10 Hz
Maximum Input:	± 10 V
Input Connection type:	miniature thermocouple
Bandwidth/Low pass filter:	-3 dB @ 10 Hz 3 pole Bessel
Secondary filters:	5, 2, 1 Hz (digital filters of e-corder)

Control Port

I²C input and output: Male and female DB-9 pin connectors.
Provides control and power.

Power requirements: ±17 V DC
+8 V DC
~3 W


Physical Configuration

Dimensions (h × w × d): 60 × 150 × 200 mm
2.4 × 5.9 × 7.9 inches

Weight: 1.5 kg (3.3 lb)

Operating conditions: 0 – 35°C
0 – 90% humidity (non-condensing)

eDAQ reserves the right to alter these specifications at any time.

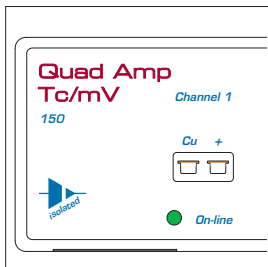

Index

B

back panel 7
block diagram 16
Bridge Amp 3

C

calibration 12
checking the Quad Tc/mV Amp 2
connecting multiple eDAQ Amps 8

E

eDAQ Amps
 Bridge Amp 3
 GP Amp 3
 Picostat 3
 Potentiostat 3
 Quad pH/mV Amp 3
 Quad Tc/mV Amp 2
 QuadStat 3

F

filtering 11
front panel 6

G

GP Amp 3

I

I²C connector 7
 pin assignments 7

L

limits of error 13

P

pH Amp
 dialog box 10
Picostat 3
Potentiostat 3
previewing the signal 10
probe calibration 12
problems 17

Q

Quad pH/mV Amp 3
Quad Tc/mV Amp 2
 setup 9–12
 use with Chart software 9
QuadStat 3

S

- setting the sensitivity 11
- signal calibration 12
- signal outputs 7
- signal preview 10
- specifications 19

T


- technical specifications 19–20
- Temperature Amp
 - dialog box 10
- thermocouple calibration 12
- thermocouple connectors 6
- thermocouple inputs 6
- thermocouple limits of error 13
- thermocouple type 11
- troubleshooting 17

U

- user modification 15

W

- warranty conditions 15, 23

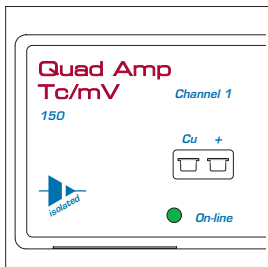

License & Warranty

Trademarks

e-corder and PowerChrom are registered trademarks of eDAQ Pty Ltd. All eDAQ software, hardware, and documentation is protected by copyright. eDAQ retains the exclusive ownership of the trademarks and registered trademarks represented by its company name, logo, and product names.

Responsibilities

You and any others using any eDAQ product agree to use it in a sensible manner for purposes for which it is suited, and agree to take responsibility for their actions and the results of their actions.

If problems arise with an eDAQ product, eDAQ will make all reasonable efforts to fix them. This service may incur a charge, depending on the nature of the problems, and is subject to the other conditions in this Agreement.

Hardware Warranty

eDAQ Pty Ltd warrants the Quad Tc/mV Amp, to be free of defects in material and workmanship for one year from the date of purchase. eDAQ Pty Ltd will repair or replace defective equipment as appropriate.

To obtain a warranty repair/replacement you must first notify us before return of the instrument and we will issue you with a RAN (return authorization number). You must ship the defective product at your

expense. We will pay return shipping. The product should be packed safely (preferably in its original packaging) and have the RAN on the shipping label. Returns sent without a RAN may be refused delivery.

This warranty does not cover hardware that has:

- been modified by the user in any way;
- been subjected to unusual physical, electrical, or environmental stress. This includes damage due to faulty power sockets, inadequate earthing, or power spikes or surges;
- been damaged because of incorrect wiring to ancillary equipment, or because of substandard, connectors or cables; or
- had the original identification marks removed or altered.

Software License

You have the non-exclusive right to use the supplied eDAQ software (Chart, Scope etc). Your employees or students, for example, are entitled to use it, provided they adhere to this agreement. Each separate purchase of the eDAQ software licenses it to be used on two computers at any given time (on one computer for data acquisition with a e-corder hardware unit, and on a second computer for the analysis of existing data files). Although multiple copies of a program may exist on several computers, more than two copies must not be used simultaneously. Departmental/company licences are

available if you wish to run more than two copies simultaneously.

Technical Support

Please register your unit to receive technical support. Technical assistance is available via email. Please describe the problem with as much detail as possible. Include a small example data file, if appropriate. Please also state:

- the model and serial number of your e-corder unit and Quad Tc/mV Amp.
- the type of computer and operating system being used (for example Windows XP, or Macintosh OS 10.2)
- the software version you are using (for example Chart v5.1.2)

We endeavor to answer all your questions, but in some cases, for example where the problem relates to the other equipment that you are using, a nominal fee may be charged.

Jurisdiction

eDAQ Pty Ltd is bound by the laws of New South Wales in Australia, and any proceedings shall be heard by the Supreme Court of New South Wales in Australia.

Disclaimer

eDAQ reserves the right to alter specifications of hardware and software without notice.

No liability can be accepted for consequential damages resulting from use of eDAQ products.